



## Pippin Pulse


# 脉冲场电泳系统



环亚生物科技有限公司 (APG BIO LTD)



美国Sage Science公司位于美国马萨诸塞州Beverly市，专注于核酸领域的样品制备。自2010年成立以来，核心产品Pippin系列全自动DNA片段回收仪因其优秀的性能表现，得到全球超过3000家客户的认可，国内装机数量400台左右。超过2万篇文献中应用，包括众多Nature、Science、Cell系列的文章已成为DNA片段筛选回收领域的行业金标准。

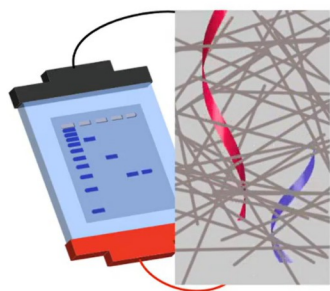


## 为什么需要脉冲场电泳？

普通琼脂糖凝胶电泳是利用分子筛的原理，分离不同大小的DNA片段。而当DNA分子大于10kb时，DNA分子的直径远远大于凝胶孔径，凝胶的分子筛就无法起到分离的作用。此时，DNA分子的迁移速率与其分子量大小不再呈现简单的相关性，因而导致无法通过观测大分子DNA的迁移速率来判断其分子量大小。

## 脉冲场电泳原理


通过快速地切换电场方向，有效地区分离大分子DNA。脉冲场电泳（PFGE, Pulsed Field Gel Electrophoresis）是一种有效分离大分子DNA的方法。通过不断改变电场方向，驱使DNA在凝胶内前向和后向来回运动，小分子DNA片段改变方向的速率要比大分子DNA片段快。

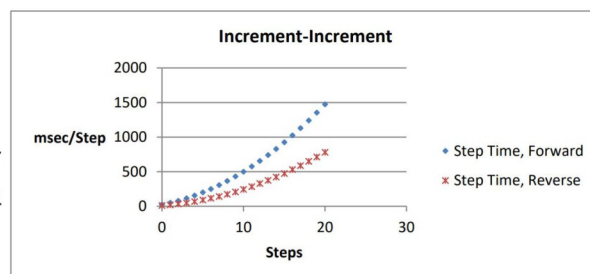

# Pippin Pulse脉冲场电泳系统

### 特点:

- ◆ 大片段DNA质控，有效分离 0.5kb-430kb;
- ◆ 预设脉冲场程序，开机即用;
- ◆ 支持自定义程序;
- ◆ 稳定耐用，加固的大直径铂电极，寿命更长
- ◆ 试剂耗材开放，成本更低

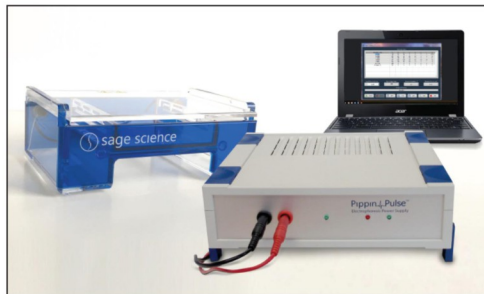

## 预设的脉冲场程序

The screenshot shows the Pippin Pulse v.1.32 application window. The 'Waveform Table' is displayed, showing various waveform types and their corresponding parameters. The table has columns for Waveform Type, V, A, B, C, D, E, F, and G. The '5-430kb' row is highlighted in blue.

| Waveform Type | V   | A   | B   | C  | D  | E  | F  | G    |
|---------------|-----|-----|-----|----|----|----|----|------|
| .5-15kb       | 100 | 90  | 36  | 0  | 0  | 0  | 0  | 5000 |
| 10-43kb       | 80  | 25  | 10  | 25 | 10 | 5  | 2  | 21   |
| 5-150kb       | 75  | 150 | 50  | 30 | 10 | 3  | 1  | 81   |
| 5-430kb       | 80  | 300 | 100 | 30 | 10 | 30 | 10 | 45   |
| 5-80kb        | 80  | 150 | 50  | 30 | 10 | 3  | 1  | 48   |
| DC Mode       | 100 | 0   | 0   | 0  | 0  | 0  | 0  | 0    |
| DC Mode 25    | 25  | 0   | 0   | 0  | 0  | 0  | 0  | 0    |

## 软件自带6种预设程序

## 支持自定义程序

| Waveform Type | V   | A  | B  | C  | D  | E | F | G    |
|---------------|-----|----|----|----|----|---|---|------|
| 0.5-15 kb     | 100 | 90 | 36 | 0  | 0  | 0 | 0 | 5000 |
| 10-48 kb      | 80  | 25 | 10 | 25 | 10 | 5 | 2 | 21   |
|               |     |    |    |    |    |   |   |      |
|               |     |    |    |    |    |   |   |      |
|               |     |    |    |    |    |   |   |      |
|               |     |    |    |    |    |   |   |      |
|               |     |    |    |    |    |   |   |      |
|               |     |    |    |    |    |   |   |      |

|                    |    |    |    |    |    |   |   |    |
|--------------------|----|----|----|----|----|---|---|----|
| 10-48 kb Version 2 | 80 | 25 | 10 | 25 | 10 | 5 | 2 | 21 |
|--------------------|----|----|----|----|----|---|---|----|

V.电压=电泳电压, 25-175 V之间。

A.运行开始时的正向时间, 1-65535毫秒。

B.运行开始时的反向时间, 1-65535毫秒。

C. 每一步增加到A的增量, 0-255毫秒。


D. 每一步增加到B的增量, 0-255毫秒。

E. 每一步增加到C的增量, 0-255毫秒。


F. 每一步增加到D的增量, 0-255毫秒。

G.每个周期的步数, 1-65535。


## 脉冲场电泳效果

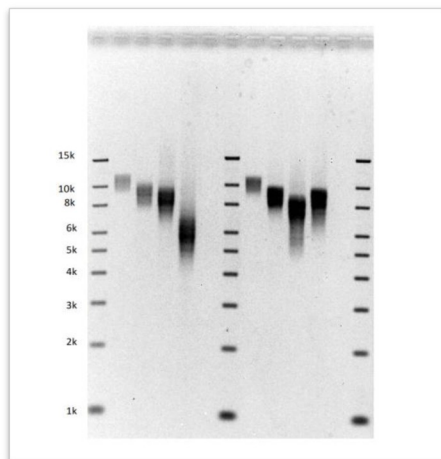


程序: 5-15 kb  
运行时间: 8 hours  
电压: 100 V  
缓冲液: 0.5×KBB  
琼脂糖胶: 0.75% Lonza SeaKem GOLD, 110ml  
Ladder: Biorad EZ Load 1kb Molecular Ruler  
# 170-8355

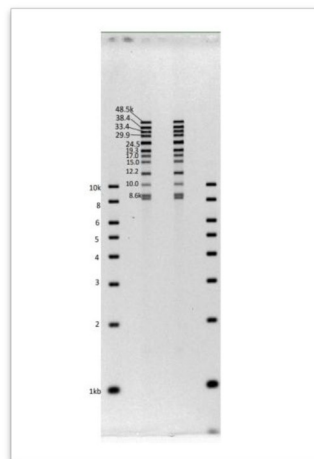


程序: 10-48 kb  
运行时间: 9 hours  
电压: 80 V  
缓冲液: 0.5×KBB  
琼脂糖胶: 0.75% Lonza SeaKem GOLD, 110ml  
Ladder: NEB Quick-load 1kb ladder #N0468S;  
Biorad CHEF DNA Size Standards 8-48 kb #170-3707

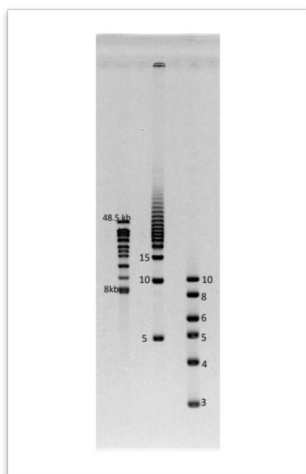

程序: 5-80 kb  
运行时间: 14 hours  
电压: 80 V  
缓冲液: 0.5×TBE  
琼脂糖胶: 1% Lonza SeaKem GOLD,  
110ml  
Ladder: Biorad CHEF DNA Size  
Standards 8-48 kb #170-3707;  
Biorad DNA Size Standard 5 Kb  
Ladder Catalog # 170-3624

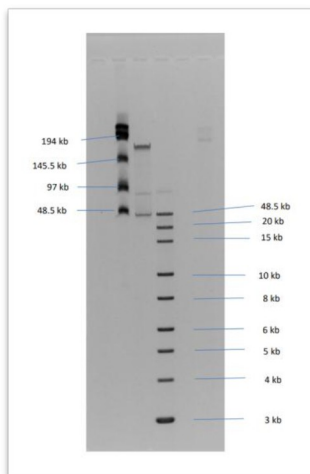

程序: 5-150 kb  
运行时间: 16 hours  
电压: 75 V  
缓冲液: 0.5×KBB  
琼脂糖胶: 0.75% Lonza SeaKem  
GOLD, 110ml  
Ladder: NEB Lambda PFG Ladder  
# N0341S; NEB 1kb Extend DNA  
Ladder #N3239

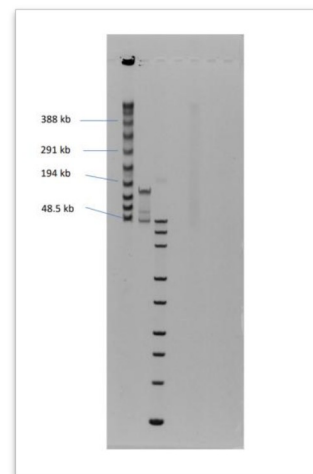

程序: 5-430 kb  
运行时间: 16 hours  
电压: 75 V  
缓冲液: 0.5×KBB  
琼脂糖胶: 0.75% Lonza SeaKem GOLD,  
110ml  
Ladder: NEB Lambda PFG Ladder #  
N0341S; Biorad CHEF DNA Size  
Standards 8-48 kb #170-3707



sage science

顶级期刊

nature

Explore content ▾ About the journal ▾ Publish with us ▾

[nature](#) > [articles](#) > [article](#)

Article | [Open Access](#) | Published: 28 April 2021

## Towards complete and error-free genome assemblies of all vertebrate species

### Isolation of high-molecular-weight DNA

**Agarose plug DNA isolation.** For tissue, HMW DNA was extracted using the **Bionano** animal tissue DNA isolation fibrous tissue protocol (cat no. RE-013-10; document number 30071), according to the manufacturer's guidelines. A total of 25–30 mg was fixed in 2% formaldehyde and homogenized using the Qiagen TissueRuptor or manual tissue disruption. For nucleated blood, 27–54  $\mu$ l was used with an adapted protocol (Bionano, personal communication) of the Bionano Prep Blood and Cell Culture DNA Isolation Kit (cat no. RE-130-10). Lysates were embedded into agarose plugs and treated with Proteinase K and RNase A. Plugs were then purified by drop dialysis with 1 $\times$  TE. DNA quality was assessed using pulse field gel electrophoresis (PFGE) (**Pippin Pulse**, **SAGE Science**, Beverly, MA) or the Femto Pulse instrument (Agilent). PFGE revealed that we isolated ultra-high-molecular-weight DNA between -100 and -500 kb long.

Nature

### Cell Reports

#### A chromosome-level reference genome and pangenome for barn swallow population genomics

##### METHOD DETAILS

**DNA extraction**  
HMW (High Molecular Weight) DNA was extracted from the muscle tissue of the samples female barn swallow with the **Bionano** animal tissue DNA isolation fibrous tissue protocol (cat no. RE-013-10; document number 30071). Approximately 55 mg of frozen muscle tissue was fixed in formaldehyde (2%) and homogenized with the Qiagen TissueRuptor. The lysate was included in agarose plugs, which were then treated with Proteinase K and RNase A. The DNA was recovered and purified from the plugs through a drop dialysis with 1 $\times$  TE. Pulsed Field Gel Electrophoresis (PFGE; **Pippin Pulse**, **SAGE Science**, Beverly, MA) and Qubit were used for DNA quality control. According to the PFGE run, a large fraction of the isolated DNA was >250Kbp.

### Cell Metabolism

#### Fatty acid overproduction by gut commensal microbiota exacerbates obesity

Graphical abstract

Authors

**Whole genome sequencing of microbes**  
E. coli TE-17 and F. intestinalis JF110941 were cultivated as described above. The culture media was centrifuged, the cell pellet was washed once with PBS, and then stored at -80°C until the DNA extraction. Bacterial genomic DNA was extracted as described above with slight modifications. Briefly, lysozyme and proteinase K were used for bacterial cell lysis. DNA quality was assessed using a spectrophotometer (DS-11, DeNovo) and Qubit fluorometer (Q33216, Thermo Fisher Scientific). For the preparation of a multiplexed microbial SMRTbell library, 1.5  $\mu$ g bacterial DNA was sheared using g-TUBE (Covaris) at 6000 rpm and purified by AMPure beads PB (Pacific Biosciences). The distribution of sheared DNA was checked by pulsed-field electrophoresis using a **Pippin Pulse** electrophoresis system (**SAGE Science**). The multiplexed microbial SMRTbell library was constructed using the "non-size selection" protocol by the manufacturer with SMRTbell Template prep kit 1.0-SPv3 (Pacific Biosciences). Equal amounts of bacterial genome libraries with barcoded adapters were pooled for sequencing. SMRTbell libraries were sequenced on the **Pacific Biosciences Sequel** platform.

Molecular Plant  
Research Article

CellPress  
OPEN ACCESS

#### The Reference Genome Sequence of *Scutellaria baicalensis* Provides Insights into the Evolution of Wogonin Biosynthesis

Qing Zhao<sup>1,2,3</sup>, Jun Yang<sup>2,3,4</sup>, Meng-Ying Cui<sup>1</sup>, Jie Liu<sup>1</sup>, Yumin Fang<sup>1</sup>, Mengxiao Yan<sup>1,2</sup>, Wenqiang Guo<sup>1</sup>, Huifen Shang<sup>1</sup>, Zhichang Xu<sup>1</sup>, Rahman Yildirim<sup>1</sup>, Jing-Xie Wang<sup>1,2</sup>, Tomáš Pluškai<sup>1</sup>, Mariette Vigouroux<sup>1</sup>, Burkhard Steuermann<sup>1</sup>, Yukun Wei<sup>1</sup>, Lei Yang<sup>1</sup>, Yonghong Hu<sup>1</sup>, Xiao-Ya Chen<sup>1,2</sup> and Cathie Martin<sup>1,2,3</sup>

Genomic DNA was extracted from leaves of a single *S. baicalensis* plant maintained in Shanghai Chenshan Botanical Garden, using a modified CTAB method (Tel-Zur et al., 1999). Quality control was done using a **SAGE Science Pippin pulse** electrophoresis system. Genomic DNA with a length of around 150 kb was sheared using a Megaruptor DNA system and the resulting fragments of 30–50 kb were collected for the

Cell子刊

nature  
biotechnology

ANALYSIS

<https://doi.org/10.1038/s41587-019-0264-0>

Corrected: Author Correction

OPEN

## Genomic and phenotypic analyses of six offspring of a genome-edited hornless bull

The integrity of the high-molecular-weight DNA samples was verified on a **Pippin Pulse gel electrophoresis system** (**SAGE Science**). The DNA was then sheared to an average size of 50 kb using a Megaruptor instrument (Diagenode) and verified on a Pippin Pulse gel. A sequencing library was prepared starting

nature  
plants

LETTERS

<https://doi.org/10.1038/s41477-019-0280-4>

## Chromosome-scale assemblies of plant genomes using nanopore long reads and optical maps

resuspended with 3 ml transposable element 10/1 buffer. The extract quality was evaluated using field inverted gel electrophoresis with the **Pippin pulse system** (**SAGE Science**). DNA samples with a fragment size above 50 kb were kept and run on BluePippin (**SAGE Science**).

nature communications

6

Article

<https://doi.org/10.1038/s41467-020-20300-3>

## A method for multiplexed full-length single-molecule sequencing of the human mitochondrial genome

2000 (Thermo Fisher Scientific) UV/Vis measurements. To determine the cell line gDNA integrity pulse-field gel electrophoresis, using the **Pippin Pulse** (**SAGE Science**) was performed. For this analysis a Sea-Kem® GOLD Agarose 1% (Lonza) gel was prepared in 0.5 $\times$  TBE buffer

Nature 子刊

SCIENCE ADVANCES | RESEARCH ARTICLE

GENETICS

## The Australian dingo is an early offshoot of modern breed dogs

Matt A. Field<sup>1,2</sup>, Sonu Yadav<sup>3</sup>, Olga Dudchenko<sup>4,5</sup>, Meera Esvaran<sup>6</sup>, Benjamin D. Rosen<sup>7</sup>, Ksenia Skvortsova<sup>8</sup>, Richard J. Edwards<sup>9</sup>, Jens Kellwagen<sup>10</sup>, Blake J. Cochran<sup>11</sup>, Bikash Manandhar<sup>9</sup>, Sonia Bustamante<sup>12</sup>, Jacob Agerbo Rasmussen<sup>13,14</sup>, Richard G. Melvin<sup>15</sup>, Barry Chernoff<sup>16</sup>, Arina Omer<sup>17</sup>, Zane Colaric<sup>18</sup>, Eva K. F. Chan<sup>2,15</sup>, Andre E. Minoch<sup>2</sup>, Timothy P. L. Smith<sup>16</sup>, M. Thomas P. Gilbert<sup>11,17</sup>, Ozren Bogdanovic<sup>2,15</sup>, Robert A. Zammit<sup>18</sup>, Torsten Thomas<sup>6</sup>, Erez L. Aiden<sup>1,5,19,20,21</sup>, J. William O. Ballard<sup>22,23\*</sup>

Copyright © 2022  
The Authors, some  
rights reserved;  
exclusive license  
American Association  
for the Advancement  
of Science. No claim  
to original U.S. Government  
Works. Distributed  
under a Creative  
Commons Attribution  
NonCommercial  
License 4.0 (CC BY-NC).

(Thermo Fisher Scientific), and molecular integrity was assessed using pulse-field gel electrophoresis. Sage Science **Pippin Pulse** assessed DNA integrity. A 0.75% KBB gel was run on the 9-hour 10- to 48-kb (80 V) program. DNA ladder used was the Invitrogen 1 Kb Extension DNA Ladder (catalog no. 10511-012). One hundred fifty nanograms of DNA was loaded on the gel.

We generated two libraries that were size-selected on Sage **BluePippin** gels (SAGE Science, Beverly, MA, USA). Libraries were sequenced on Sequel machines with 2.0 chemistry recording 10-hour movies. Sequencing was conducted at the Arizona Genomics Institute, University of Arizona.

Science Advances

PACBIO®

## Procedure & Checklist - Using the Sage Science™ Pippin Pulse Electrophoresis Power Supply System

PacBio官方推荐Pippin Pulse

## 代表文献

### Pippin Pulse + Nanopore

- Long, Meng, et al. "Comparative genomic analysis provides insights into taxonomy and temperature adaption of *Aeromonas salmonicida*." **Journal of Fish Diseases** 46.5 (2023): 545-561.
- Keraite, Ieva, et al. "A method for multiplexed full-length single-molecule sequencing of the human mitochondrial genome." **Nature Communications** 13.1 (2022): 5902.
- Ong, Chian Teng, et al. "Adaptive sampling during sequencing reveals the origins of the bovine reproductive tract microbiome across reproductive stages and sexes." **Scientific reports** 12.1 (2022): 15075.
- Wunderlich, Stephanie, et al. "Targeted biallelic integration of an inducible Caspase 9 suicide gene in iPSCs for safer therapies." **Molecular Therapy-Methods & Clinical Development** 26 (2022): 84-94.
- Russo, Alessia, et al. "Low-input high-molecular-weight DNA extraction for long-read sequencing from plants of diverse families." **Frontiers in Plant Science** 13 (2022): 1494.
- Player, Robert A., et al. "A novel *Canis lupus familiaris* reference genome improves variant resolution for use in breed-specific GWAS." **Life science alliance** 4.4 (2021).
- Young, Amy E., et al. "Genomic and phenotypic analyses of six offspring of a genome-edited hornless bull." **Nature biotechnology** 38.2 (2020): 225-232.
- Belser, Caroline, et al. "Chromosome-scale assemblies of plant genomes using nanopore long reads and optical maps." **Nature plants** 4.11 (2018): 879-887.

### Pippin Pulse + PacBio

- Takeuchi, Tadashi, et al. "Fatty acid overproduction by gut commensal microbiota exacerbates obesity." **Cell Metabolism** (2023).
- Hadzhiev, Yavor, et al. "The miR-430 locus with extreme promoter density forms a transcription body during the minor wave of zygotic genome activation." **Developmental Cell** 58.2 (2023): 155-170.
- Kurokuchi, Hiroyuki, et al. "Telomere-to-telomere genome assembly of matsutake (*Tricholoma matsutake*)." **DNA Research** 30.3 (2023): dsad006.
- Hixson, Kim K., et al. "Annotated genome sequence of a fast-growing diploid clone of red alder (*Alnus rubra* Bong.)." **G3: Genes, Genomes, Genetics** (2023): jkad060.
- Field, Matt A., et al. "The Australian dingo is an early offshoot of modern breed dogs." **Science Advances** (2022): eabm5944.
- Pootakham, Wirulda, et al. "A Chromosome-Scale Genome Assembly of *Mitragyna speciosa* (Kratom) and the Assessment of Its Genetic Diversity in Thailand." **Biology** 11.10 (2022): 1492.
- Rahman, M. Julhasur, et al. "LINE-1 retrotransposons facilitate horizontal gene transfer into poxviruses." **Elife** 11 (2022): e63327.
- Zhao, Qing, et al. "The reference genome sequence of *Scutellaria baicalensis* provides insights into the evolution of wogonin biosynthesis." **Molecular plant** 12.7 (2019): 935-950.

### Pippin Pulse + Bionano

- Secomandi, Simona, et al. "A chromosome-level reference genome and pangenome for barn swallow population genomics." **Cell Reports** 42.1 (2023): 111992.
- Dahn, Hollis A., et al. "Benchmarking ultra-high molecular weight DNA preservation methods for long-read and long-range sequencing." **GigaScience** 11 (2022).
- Palmada-Flores M, Orkin J D, Haase B, et al. A high-quality, long-read genome assembly of the endangered ring-tailed lemur (*Lemur catta*) [J]. **GigaScience**, 2022, 11.
- Canaguier A, Guilbaud R, Denis E, et al. Oxford Nanopore and Bionano Genomics technologies evaluation for plant structural variation detection [J]. **BMC genomics**, 2022, 23(1): 1-17.
- Shukla H, Suryamohan K, Khan A, et al. Near-chromosomal de novo assembly of Bengal tiger genome reveals genetic hallmarks of apex-predation [J]. **bioRxiv**, 2022.
- Rhie, Arang, et al. "Towards complete and error-free genome assemblies of all vertebrate species." **Nature** 592.7856 (2021): 737-746.
- Belser, Caroline, et al. "Telomere-to-telomere gapless chromosomes of banana using nanopore sequencing." **Communications biology** 4.1 (2021): 1047.
- Zheng, Yuting, et al. "Rapid prenatal diagnosis of Facioscapulohumeral Muscular Dystrophy 1 by combined Bionano optical mapping and karyomapping." **Prenatal Diagnosis** 40.3 (2020): 317-323.

## 参考客户

华大基因  
BGI

Novogene  
诺禾致源

ANOROAD  
安诺优达

组 NextOmics 未来组  
组 GrandOmics 希望组

中国科学院昆明动物研究所  
KUNMING INSTITUTE OF ZOOLOGY, CAS

复旦大学

中山大学  
SUN YAT-SEN UNIVERSITY

中国医学科学院北京协和医学院  
Chinese Academy of Medical Sciences & Peking Union Medical College

北京大学

天津大学  
TIANJIN UNIVERSITY

军事医学科学院  
ACADEMY OF MILITARY MEDICAL SCIENCES

中国农业大学  
China Agricultural University

首都医科大学  
CAPITAL MEDICAL UNIVERSITY

## Sage Science片段回收系列



### Pippin Prep™

常规通用型

回收范围: 100bp - 1.5kb  
样本数量: 5个



### BluePippin™

全长型

回收范围: 100bp - 50kb  
样本数量: 5个



### PippinHT™

高速高通量型

回收范围: 100bp - 50kb  
样本数量: 24个



### Pippin 24™

多功能型

片段筛选回收、电泳图谱分析、  
核酸浓度测定等  
样本数量: 24个



### SageELF™

多通道回收型

回收范围: 100bp - 40kb  
样本数量: 2个



### HLS2

超大片段型

回收范围: 高达2Mb  
样本数量: 4个



## 环亚生物科技有限公司

美国Sage Science大中华区独家代理商



扫码关注微信公众号

地址: 上海市松江区莘砖公路518号9幢502室

电话: 021-54583565

邮箱: [info@apgbio.com](mailto:info@apgbio.com)

网站: [www.apgbio.com](http://www.apgbio.com)